UYA 4Y SERIES MICROBALANCES

release date: 01-07-2016









APPLICATION

4Y series is a modern weighing device, especially useful when the measurement requires perfect accuracy and high speed.

FEATURES OF 4Y SERIES MICROBALANCE









UYA 2.4Y (standard solutions for any operating conditions)






UYA 2.4Y (control of piston pipettes volume)

ACCURACY

Measurement accuracy is guaranteed with an adjustment procedure carried out using an internal adjustment weight. This fully automatic process is controlled by a module intended to diagnose ambient conditions change (on-line). Adjustment processes (internal and external) can be performed in accordance with a specially designed schedule.

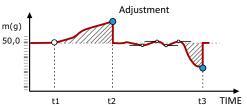

FUNCTIONALITY

4Y series microbalance is an optimized modern device which features an option of automatic level control (Level SENSING) as one of numerous functions.

The practical effect of balance customization are individual user profiles and gradable permission levels for access to balance menu.

Programmable proximity sensors offer wide range of possibilities: weighing chamber control, zeroing, tarring, printout.

Several functions such as differential weighing facilitate multi-stage mass control of one and the same sample subjected to various processes. Pipettes calibration function is an ergonomic tool designed to calibrate and control piston pipettes with the use of gravimetric method. MEDIA module, as one of the greatest 4Y series assets, provides the user with an on-screen help and support.

COMPLIANCE WITH REGULATIONS

Owing to security system and possibility to document the process by means of printouts (standard/editable), the 4Y balance meets requirements imposed by GLP/GMP systems for various industries (pharmacy, petrochemistry, environmental protection).



PARTS COUNTING



DOSING



CHECKWEIGHING



FORMULATION



PERCENT WEIGHING [%]



STATISTICS



ANIMAL WEIGHING



DIFFERENTIAL WEIGHING



PIPETTES CALIBRATION



SQC



AUTOTEST



AUTOMATIC DOOR



AIR DENSITY CORRECTION



GLP / GMP PROCEDURES



INFRARED SENSORS



AMBIENT COND. MONITORING



UNITS

MEDIA module support within a reach of your hand

A sheer novelty of 4Y microbalance is the MEDIA module.
With it you can learn using videos providing information on any relevant matter.




MEDIA module is a convenient way allowing you to broaden your knowledge on mass measurement but noe exlusively (recommendations, SOP, reminders, self-designed testing procedures).

	UYA 2.4Y	UYA 2.4Y.F
	M 16	M 16
Max load	2,1 g	2,1 g
Min load	0,01 mg	0,01 mg
Readability	0,1 µg	0,1 µg
Repeatability	0,3 μg (Rt≤0,2g)	0,4 µg (Rt≤0,2g)
	0,6 µg (0,2g <rt≤2g)< td=""><td>0,6 µg (0,2g<rt≤2g)< td=""></rt≤2g)<></td></rt≤2g)<>	0,6 µg (0,2g <rt≤2g)< td=""></rt≤2g)<>
Linearity	±1,5 μg	±1,5 μg
Eccentric load deviation	1,5 µg	1,5 µg
Sensitivity offset	1,5 × 10 ⁻⁶ × Rt	1,5 × 10 ⁻⁶ × Rt
Sensitivity temperature drift	1 × 10 ⁻⁶ / °C × Rt	1 × 10 ⁻⁶ / °C × Rt
Sensitivitytime drift	1 × 10 ⁻⁶ / Year × Rt	1 × 10 ⁻⁶ / Year × Rt
Minimum weight (USP)	0,8 mg	0,8 mg
Minimum weight (U = 1%, k = 2)	0,08 mg	0,08 mg
Pan size	ø 16 mm	ø 50 mm
Weighing chamber dimensions	ø 90 × 90 mm	ø 118 × 35 mm
Stabilization time	10÷20 s	
Adjustment / Calibration	automatic (internal)	
Power supply	13,5 ÷ 16 V DC / 2,1 A	
Casing of the terminal	ABS plastic	
Display	colour 5,7"(640x480) with a resistive touch screen	
Processor	2 × 1 GHz	
Memory	RAM: 256 MB DDR2, flash: 8 GB microSD	
Interface	2×USB host, 2×RS 232, Ethernet 10/100 Mbit, Wireless Module - optional	
Audio module	YES (voice messages support)	
Video support	YES (videos and multimedia instructions)	
IN / OUT	4 in / 4 out (digital)	
Ambient conditions:		
Working temperature	+10 ° ÷ +30 °C	
Change rate of working temperatures	±0,3°C/1h (±0,5°C / 12h)	
Relative air humidity **	40% ÷ 80%	
Change rate of atmospheric humidity	40% ÷ 60% (±5% / 4h)	

Rt - net weight

Dimensions:

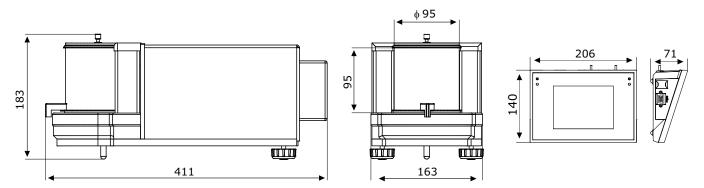

Owing to fastidious selection of mechanical design components our balance allows to obtain great stability and repeatability of indications regardless of measured sample size and conditions of use. Multi-shield mechanical design of the weighing module provides excellent thermal stability. Even with challenging ambient temperature your weighing is precise and accurate.


4Y microbalances offer modern hardware and software. In-build programs, Windows Embedde Compact 7 operating system, Flash memory, Double Hardware system guarantee fast measurement and reliability when it comes to data acquiring and


processing. The 4Y series comprises 5,7" colour touchscreen providing ever more increased balance operation functionality and even more practical results presentation. Complex databases allow measurement record along with printout and export option.


^{** -} Non-condensing conditions

Dimensions:

F version

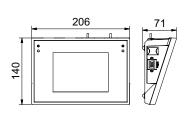

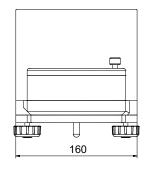

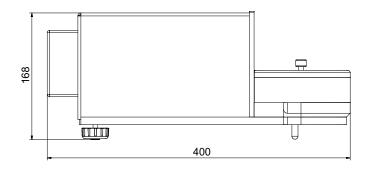

Anti vibration table for microbalance	Ambient conditions module	
Profesional weighing table	Additional LCD display "WD-5"	
Impact Epson printer	PC keyboard	
Label printer Citizen	Power adapter with battery and charger ZR-02	
Anti draft shield for microbalances	Mass standard	
"Tare" and "Print" foot button	Antistatic cable	
"PW-WIN" computer software	Bar code scanner	
"RAD-KEY" computer software	Cable RS 232 (balance - computer) "P0108"	
"REC-FS" computer software	Cable RS 232 (balance - Epson, Citizen printer) "P0151"	
Antistatic ionizer DJ-03		